

updater.zip

Contains files used to update the device and the PC software.

Procedure:

1) Unzip the zip file. You get a folder "updater" and a program file "trace_mon.exe". The updater is designed to update the unit and the program file is the respective software for the control PC.

2) Copy the folder "updater" in the root directory of a USB stick

3). Insert the USB stick into the BioPAT ® Trace and turn on the device. After booting (white LED), the LED changes to red. Turn it off again and repeat the process until after booting a traffic light signal red, yellow, green appears.

(Note: The traffic signal serves as verification that the USB stick has been detected properly. Sometimes USB memory sticks boot too slowly and it may happen that the BioPAT ® Trace does not realize it has a USB Stick with further instructions on board. In this case several attempts are necessary.)

4) To update the program file overwrite the existing file "trace_mon.exe" with the new version in your program directory on your PC.